



863 Valley View Road, Eighty Four PA 15330
724-941-9701 • skcinc.com

BioSampler Case with Mounting Rod Operating Instructions

Assembly

1. Remove nylon harness, clip (with thumbscrew), and mounting rod from case (mounting rod is between back of foam insert and back of case).
2. Slide large end of mounting rod through the larger loop on the nylon harness until the harness loop rests in the indentation on the mounting rod.
3. Slide the small end of the mounting rod into the small hole on the clip.
4. Tighten the thumbscrew.
5. Screw threaded end of mounting rod into the semi-circular hub on top of the case.
6. Place BioSampler (assembled) into the indentation in the hub.
7. Loosen the thumbscrew on the clip and place clip indentation around the BioSampler nozzle. Tighten the thumbscrew.
8. Wrap nylon harness around the BioSampler and secure Velcro fastener.

Disassembly

1. Remove BioSampler and unscrew mounting rod.
2. Remove nylon harness.
3. Loosen the thumbscrew and remove clip.
3. Place nylon harness and clip in the back, right foam compartment.
4. Place the mounting rod between the back of the foam insert and the back of the case.

Caution: Failure to replace parts in the proper foam compartments will cause distortion to the inside of the carry case and reduce its ability to protect glass components.