



### Surface SWYPE for Aromatic Amines Cat. No. 769-1021

1. Lightly spray the area or item (e.g., workbench, tool, control knob) with Cleaning/Developing Solution for Aromatic Amines Cat. No. 769-1041 or 769-1061.
2. Wait at least 30 seconds and then wipe with a Surface SWYPE pad.
3. Allow 3 minutes for the color to develop on the pad. Aromatic amines will produce the following colors:
  - *MOCA and MDA*: Red-orange
  - *Ethacure*: Blue
  - *Phenylene diamine*: Purple

**Notes:** *Although the color is stable, it is recommended that you read the color after 3 minutes.  
Known interference: MDI*

*To request testing of other specific aromatic amines, contact SKC.*

#### **SKC Limited Warranty and Return Policy**

SKC products are subject to the SKC Limited Warranty and Return Policy, which provides SKC's sole liability and the buyer's exclusive remedy. To view the complete SKC Limited Warranty and Return Policy, go to [skcinc.com/warranty](http://skcinc.com/warranty).